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## Experiment 8 Qualitative determination of Carbohydrates

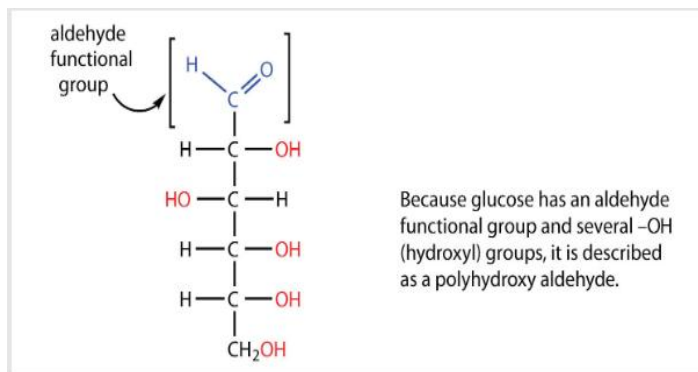
This experiment is intended to introduce you to one of the three major classes of macronutrients found in food, carbohydrates. To identify carbohydrates (sugars) as being polyhydroxylated aldehydes and ketones. Mono or di- or polysaccharides and their chemical characteristics. You will also learn a variety of ways to categorize carbohydrates and several tests used in the analysis of carbohydrates.

### Introduction of Carbohydrate

**Carbohydrate** is the primary source of energy for brain, retina and erythrocytes and so the glucose concentration requires precise regulation that is accomplished by many metabolic processes and interrelated with several glands and organs. If any disturbance occurs to the glands or organs, a loss of glycemic control will occur which results in the formation of a pathological state. However, the **diabetes mellitus** is the result of glucose metabolism abnormality, and is often has different causes, which include genetic, autoimmune and other causes.


All carbohydrates consist of carbon, hydrogen, and oxygen atoms and are polyhydroxy aldehydes or polyhydroxyketones or are compounds that can be broken down to form such compounds. Examples of carbohydrates include glucose, ribose, fructose (fruit sugar), sucrose (table sugar), lactose (milk sugar), maltose (malt sugar), galactose, starch, and structural materials such as cellulose. The molecular formula for a carbohydrate can be considered as  $C_n(H_2O)_n$ , where n is the number of carbon atoms. Example, because its formula is  $C_6H_{12}O_6$ , glucose was once thought to be a "carbon hydrate" with the structure  $C_6(H_2O)_6$ .

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**Fructose has a ketone group**, so it is considered a **polyhydroxy ketone**.

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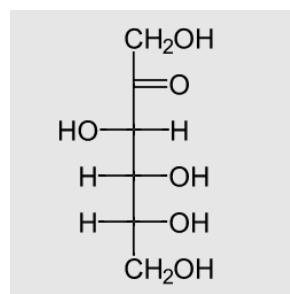## Classification according to Molecular size or Complexity:

- A- Simple Carbohydrates.** often referred to as **simple sugars or monosaccharides**, contain one unit of saccharide and cannot be broken down into smaller carbohydrates, such as glucose, fructose, galactose, and mannose.

Monosaccharide can be also classified according to different factors as follow:

### a. Classification according to **Number of carbon atoms**

1. **Triose C3 sugars**, example: dihydroxyacetone and glyceraldehydes.
2. **Tetrose C4 sugars**, example: D-erythrose, D-threose, and D-erythrulose.
3. **Pentose C5 sugars**, example: nucleic acids and coenzymes NAD, Ribose that is a structural element of ATP and flavor proteins: Arabinose, ribulose, and xylose.
4. **Hexose C6 sugars**, example: glucose, fructose.

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### b. Classification according to **Functional group**

1. **Aldose**: sugars having an aldehyde functional group R-HC=O, example: glucose
2. **Ketose**: sugars having a ketone functional group R<sub>2</sub>-C=O, example: fructose.

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### c. Classification according to **Reactivity in Redox Reactions**

1. **Reducing sugars** oxidized by Tollens' reagent or Benedict's or Fehling's reagents). Examples are glucose, fructose, glyceraldehydes, lactose, arabinose and maltose
2. **Non-reducing sugars** not oxidized by Tollens' reagent or Benedict's or Fehling's reagents. Example: sucrose, starch.

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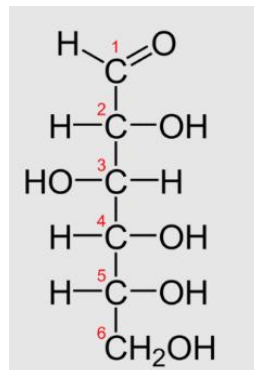
for example, glucose is an aldose (it contains an aldehyde functional group) and glucose is also a hexose (it contains six carbon atoms). Therefore, glucose is considered as an aldohexose.

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Glucose is classified as an aldohexose, because it contains six carbon atoms and an aldehyde functional group.

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**Complex Carbohydrates.** Those containing more than one group of saccharides are complex carbohydrates. Through a process known as hydrolysis, complex carbohydrates can be broken down into smaller carbohydrate units.

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1. **Disaccharides** contain two units of monosaccharide., Hydrolysis can break disaccharides into two monosaccharide units and lose one molecule of water. such as:

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- **Maltose:** two glucose residues in 1, 4 linkages.

- **Lactose** can be found in milk. On hydrolysis, it produces D-galactose and D-glucose. it is a reducing disaccharide

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- **Sucrose:** Cane sugar (table sugar) or sucrose is a disaccharide of fructose and glucose.

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2. **Oligosaccharides** have 3-6 units of monosaccharides, Oligosaccharides can be broken by hydrolysis into 3-6 monosaccharide units and lose one or more water molecules.

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3. **Polysaccharides** (hundreds or thousands of units), Polysaccharides can be broken into 7 or more units of monosaccharides by hydrolysis and one or more molecules of water can be lost. Such as:

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-**Starch** is a storage form of energy in plants. It contains two polymers composed of glucose units: amylose (linear) and amylopectin (branched). The complete hydrolysis of starch yields, in successive stages, glucose:

starch → dextrins → maltose → glucose

In the human body, several enzymes known collectively as amylases degrade starch sequentially into usable glucose units.

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-**Glycogen** is a storage form of energy in human and animals. It is a branched polymer composed of glucose units. It is more highly branched than amylopectin. glycogen is found as granules in liver and muscle cells. When fasting, animals draw on these glycogen reserves during the first day without food to obtain the glucose needed to maintain metabolic balance.

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-**Cellulose** is a structural polymer of glucose units found in plants. It is a linear polymer with the glucose units linked through β-1,4-glycosidic bonds.

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## Qualitative tests for carbohydrates:

There are several difficulties in their qualitative as well as quantitative analysis when analyzing a sample containing a mixture of carbohydrates, particularly sugar. These problems are attributed to their structural and chemical resemblance, as well as their stereoisomerism. It is therefore necessary to determine, during biochemical investigations, whether or not a given sample contains carbohydrates. In order to establish the presence or absence of a sugar or a carbohydrate in a sample, several rapid tests are available. Such tests are based on specific typical color reactions for their group. The sensitivity of these tests can be confirmed by using different concentrations of sugar solutions (0.1- 1 percent ).

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For all carbohydrates, whether free or in combined form, it is a group test. It is routinely used to detect the presence of carbohydrates, despite its limitations.

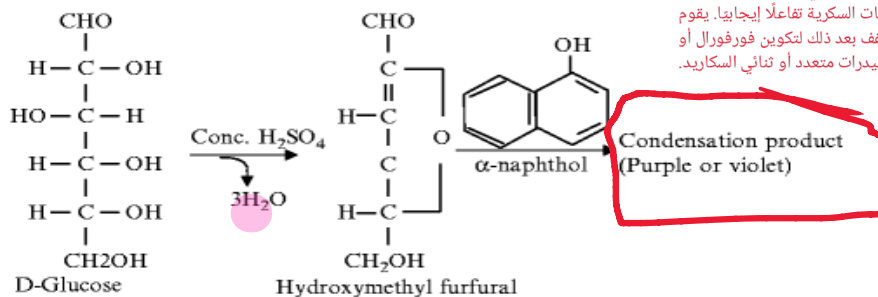
In order to detect the presence of carbohydrates in a solution, the most common tests used are:

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The response is based on the fact that H<sub>2</sub>SO<sub>4</sub> concentrated catalyzes the dehydration of sugars to form furfural (from pentoses) or hydroxymethyl furfural (from hexoses). These furfurals then condense to give a purple or violet colored product with sulfonated alpha-naphthol. A+ve reaction is also provided by polysaccharides and glycoproteins. The acid first hydrolyses it into monosaccharides, which are then dehydrated to form furfural or its derivatives, in the case of the carbohydrate being a poly or disaccharide.

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## B. Differentiation test of Carbohydrates

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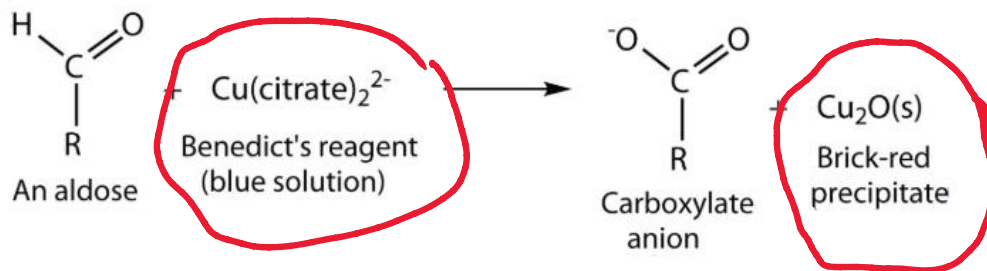
Benedict's test determines whether a **monosaccharide or disaccharide is a reducing sugar**. To give a **positive test**, the carbohydrate must **contain either a free aldehyde group or a ketone group**. Benedict's reagent is an **alkaline solution containing cupric ions**, which **oxidize the aldehyde to a carboxylic acid**. In turn, **the cupric ions are reduced to cuprous oxide**, which forms a **red precipitate**.

Give **positive result with all reducing sugar as glucose, fructose, lactose, maltose, galactose, except for sucrose and starch.**

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Consist of **cupric ion in acidic solution (acetic acid)**. This test is performed to differentiate between **mono- and disaccharide reduction**. Monosaccharides are more reactive reducing agents than disaccharides and thus react in about 1-2 minutes, while it takes 7-12 minutes for the reducing disaccharides to get hydrolysed and then react in the acidic solution. Therefore, it is possible to detect the difference in property reductions.

It gives positive red precipitate result with glucose and fructose and other monosaccharides.

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This test is a **keto hexose-specific timed colour reaction**. It is therefore used for the differentiation of **aldoses from ketoses**. Dehydration is carried out in the presence of **concentrated HCl or H2SO4 with resorcinol**, **keto hexoses to yield 4-hydroxy methyl furfural** more rapidly than aldohexoses. In addition, these furfural derivatives condense to form a **red coloured complex with resorcinol**.

It gives positive result with fructose. And in **prolong heating it will give a positive result with sucrose also due to hydrolysis to fructose and glucose, while other sugar gives a negative result.**

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### 4. Iodine test:

Iodine forms color complex with starch, glycogen and dextrins. The color varies according to the **chain length and the size of the molecule**. **Starch gives blue color, glycogen gives purple** while **dextrins forms red to colorless complexes**. The complexes with iodine are very unstable. The iodine is readily removed from these compounds by alcohol, NaOH, Na2S2O3 or heating. This indicates the **weak interaction between iodine and these compounds**.

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### 1. Molisch's test:

#### Procedure

1. In a test tube, add 2 ml of test solution ( unknown)
2. Add 2 drops of Molisch's reagent and mix by gentle shaking
3. Tilt the tube and carefully pour 1 ml conc. H<sub>2</sub>SO<sub>4</sub> along the side of the tube (no mixing)
4. Violet ring at the junction of the two liquids indicates a positive result.

*Note: The appearance of purple or violet ring confirms the presence of carbohydrate.*

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**Interpretation:** The formation at the junction of two layers of a **purple or violet ring or zone** indicates the presence of **carbohydrates**.

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#### Precautions:

- (i) The solution of alpha-naphthol is unstable and should be made fresh.
- (ii) Conc.H<sub>2</sub>SO<sub>4</sub> should be carefully added along the sides of the test tube, causing the contents of the tube to be minimally disturbed.

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### 2. Iodine test:

1. To two ml of your test solution ( unknown), add few drops of diluted HCl
2. Add 2 drops of iodine solution
3. A **blue color** indicates positive result
4. Heat the colored solution and observe the disappearance of the color

Iodine test give A+e result ( blue color complex) with starch as a polysaccharides.

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### 3. Benedict's test

1. In a tube add 5 ml of benedict reagent
2. Add 8 drops of your test solution (unknown) and mix
3. Heat the tube in boiling water bath for 5 min
4. Positive result is indicated by the **change in color or the formation of a precipitate yellow to brick red color**, depending upon the conc. of the glucose or reduced sugar in the solution.

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*Note: The appearance of red precipitate confirms the presence of carbohydrates.*

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### 4. Barfoed's test

1. Add 1 mL of the test solution (unknown) to a test tube.
2. Add 5 ml Barfoed's reagents to test tube and mix
3. Place the test tubes in a water bath and observe the time when a precipitate appears.

Thin red precipitates indicate the presence of a **reduction of monosaccharide** at the bottom or sides of the tube.

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**Positive Barfoed's test for monosaccharides: glucose, fructose.**

**Note:** The boiling should not be prolonged beyond 1-2min, otherwise the disaccharide reduction will respond to this test as well.

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## 5. Seliwanoff's test

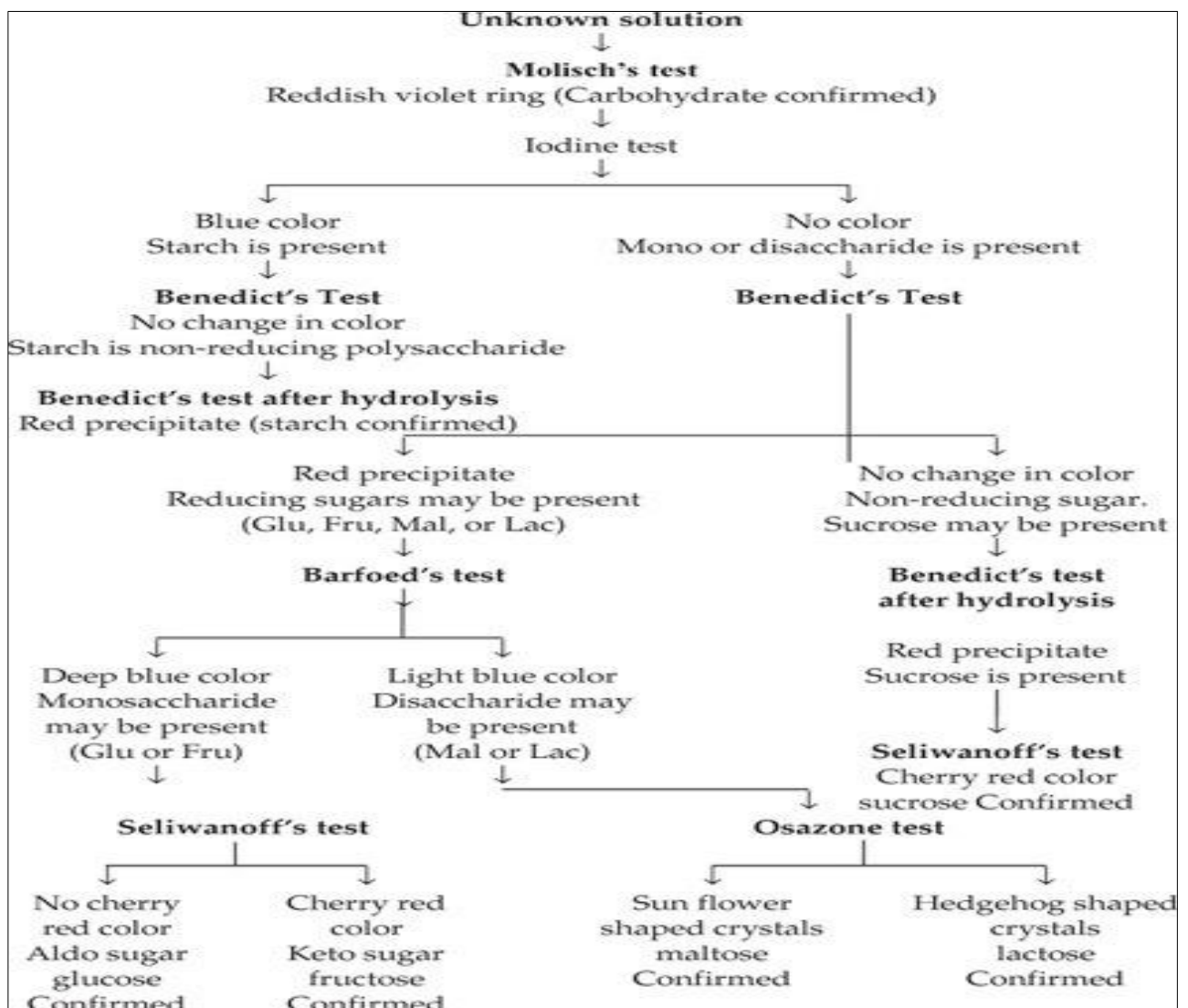
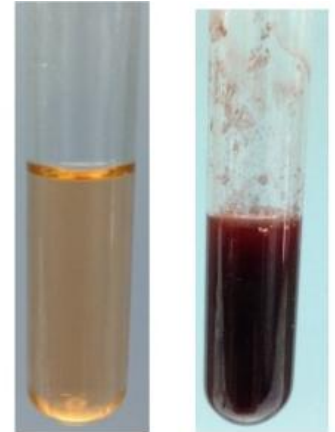
1. Add 1 mL of the test solution (unknown) to about 2 mL of the Seliwanoff reagent
2. Heat the test tube in boiling water bath for 10-20 min or just heat to boiling.

A positive reaction is indicated by a **deep red color and formation of brown precipitate**.

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(The presence of **ketohexose** indicates the appearance of a **red colour (fructose)**).

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## Summary

Biochemical test	Glucose	Fructose	Maltose	Lactose	Sucrose	Starch
Molisch	<b>Positive</b>	<b>Positive</b>	<b>Positive</b>	<b>Positive</b>	<b>Positive</b>	<b>Positive</b>
Benedict	<b>Positive</b>	<b>Positive</b>	<b>Positive</b>	<b>Positive</b>	<b>Negative</b>	<b>Negative</b>
Barfoed	<b>Positive</b>	<b>Positive</b>	<b>Negative</b>	<b>Negative</b>	<b>Negative</b>	<b>Negative</b>
Seliwanoff	<b>Negative</b>	<b>Positive</b>	<b>Negative</b>	<b>Negative</b>	<b>Positive</b>	<b>Negative</b>
Iodine	<b>Negative</b>	<b>Negative</b>	<b>Negative</b>	<b>Negative</b>	<b>Negative</b>	<b>Positive</b>